Existent Gum in Fuels by Jet Evaporation

test method

Gum formed during fuel storage can deposit on induction system surfaces, intake valves, stems and guides. To test for gum content, a 50mL sample is evaporated in an aluminum block bath for a specified period under controlled conditions of temperature and flow of air (aviation and motor gasolines) or steam (aircraft turbine fuel).

existent gum test apparatus

Evaporates aircraft turbine fuel and motor and aviation gasoline samples under controlled conditions in accordance with ASTM specifications. Consists of a high temperature evaporation bath with 100mL test beakers and, for aircraft turbine fuels, a steam generator and steam superheater.

evaporation baths

- · Conforming to ASTM D381 and related specifications
- Choice of three-unit and six-unit models
- Safety top assembly allows for easy positioning and connection of all conical adapters to the bath
- · Digital flowmeter for accurate and precise air flow measurement
- Built-in steam superheater
- Microprocessor programmable high accuracy temperature control
 Built-in pressure regulator

Electrically heated baths for determining existent gum in aircraft turbine fuels by steam-jet evaporation and in motor and aviation gasolines by air-jet evaporation. Fully insulated, aluminum block design assures safe, efficient high temperature operation. Equipped with air/steam pressure regulator with gauge and a digital flowmeter for adjusting air flow per ASTM specifications. Stainless steel jets deliver air or steam flow to the test wells through removable brass conical adapters. Microprocessor PID control provides quick temperature stabilization without overshoot, and the bath is protected by an overtemperature control circuit that interrupts power should bath temperature exceed a programmed cut-off point. Dual LED displays provide actual and setpoint temperature values in $^{\circ}C/^{\circ}F$ format. Communications software (RS232, etc.), ramp-to-set and other enhanced features are available as extra cost options. Contact your Koehler representative for information.

Model **K33900** with Built-in Superheater–Six-unit bath with a builtin thermostatically controlled superheater which delivers dried steam to the bath inlet for steam-jet method testing of aircraft turbine fuels. Has digital indicating solid state bath temperature control with digital setpoint and display.

Model **K33780/K33781** – Three-unit bath without built-in superheater. All controls are housed in the bath cabinet.

ordering information

catalog no. description

- K33900 Existent Gum Evaporation Bath, 6-Unit with Superheater, 220-240V 50/60Hz
 K33780 Existent Gum Evaporation Bath, 3-Unit, 115V 60Hz
- **K33781** Existent Gum Evaporation Bath, 3-Unit, 220-240V 50/60Hz



K33900 Existent Gum Evaporation Bath

specifications

Conforms to the specifications of: ASTM D381; IP 131; IP 540; ISO 6246; DIN 51784; FTM 791-3302; NF M 07-004

Testing Capacity

K33900: 6 sample beakers K33780 and K33781: 3 sample beakers Maximum Temperature: 475°F (246°C) Temperature Control Stability: 1°F (0.5°C) Bath Configuration: machined aluminum block with multiple cartridge heaters

Heater Range

K33900: 0-3000W K33780 and K33781: 0-1500W Superheater: (Model K33900 only) Superheating chamber and condensate trap constructed of stainless steel. Solid state thermoregulator (0-550°F) Heater Range: 0-3000W

Electrical Requirements

K33900: 220-240V 50/60Hz, Single Phase, 30A K33780: 115V 60Hz, Single Phase, 13.0A K33781: 220-240V 50/60Hz, Single Phase, 6.8A

Included Accessories Conical Brass Adapters for air/steam jets

Dimensions Ixwxh,in.(cm) K33900 21"x20.5"x28" (53.34x52.07x71.12) Net Weight 230 lbs (104.3kg) K33780/K33781 32.5x11x19 (83x28x48) Net Weight 85 lbs (38.6kg) Shipping Information: Shipping Weight/Dimension K33900 313 lbs (142kg) / 17.2 Cu. ft. K33780/K33781 140 lbs (63.5kg) / 8.3 Cu. ft.



85 Corporate Drive, Holtsville, New York 11742 1-800-878-9070 (in u.s. only) TEL: +1 631 589 3800 FAX: +1 631 589 3815 Email: sales@koehlerinstrument.com www.koehlerinstrument.com

Existent Gum in Fuels by Jet Evaporation

steam generator

- · For steam-jet method testing of aircraft turbine fuels
- Meets output requirements of three-unit and six-unit evaporation baths
- Electrically heated for clean, efficient operation and ease of installation

•Meets applicable ASME, NEC standards; UL listed, CSA approved Electrically heated boiler provides instantaneous and reserve steam capacity for steam-jet evaporation tests. Easy to install and operate; electrical heating eliminates the need for on-site fuel combustion. Requires only a water feed source and electrical hook-up. Ruggedly constructed, with long life industrial grade incoloy heating element. Includes a full range of safety features: automatic water level control and low water cut-off; steam safety valve; high-limit pressure cut-out with manual reset; steam pressure gauge.

specifications

Output: 54.1 lbs steam/hr at 212°F Bhp Rating: 1.83 kW Rating: 18

Dimensions lxwxh,in.(cm) 20x28x36 (51x71x91) Net Weight: 185 lbs (83.9kg)

Shipping Information

K33850/480603

Shipping Weight: 200 lbs (91kg) Dimensions: 18 Cu. ft.

ordering information

catalog no. description K33850 Steam Boiler, 120/240V 60Hz, Three Phase K33850/208601 Steam Boiler, 208V 60Hz, Single Phase, 87A K33850/208603 Steam Boiler, 208V 60Hz, Three Phase, 50A K33850/240601 Steam Boiler, 240V 60Hz, Single Phase, 75A K33850/240603 Steam Boiler, 240V 60Hz, Three Phase, 43A Steam Boiler, 380V 50/60Hz, Three Phase, 27A K33850/380603 K33850/415503 Steam Boiler, 415V 50Hz, Three Phase, 25A

Other electrical configurations for the Steam Boiler are available. Please inquire with Koehler Customer Service for additional information.

Steam Boiler, 480V 60Hz, Three Phase, 22A



accessories

catalog no.	description	qty	
K33710	Sample Beaker, 100mL spun copper, 50x78mm	6	
332-002-017	Sample Beaker, Borosilicate Glass, 100mL		
250-000-03F	ASTM 3F Thermometer Range: 20 to 760°F	2	
250-000-03C	ASTM 3C Thermometer Range: -5 to +400°C		
K33810	Steam Superheater		
	Provides dry superheated steam for evaporation		
	baths not equipped with a built-in superheater. Use		
	together with an outside steam source for steam-jet		
	method testing of aircraft turbine fuels. Superheating		
	chamber and condensate trap are constructed entirely of		
	stainless steel. Solid state temperature controller adjusts		
	between 0-550°F. Equipped with steam inlet and outlet		
	connections and condensate drain valve. Steel exte	erior	
	has a chemical resistant polyurethane enamel finish		

Dimensions 5x27x9.5" (13x70x24cm). Shipping Weight: 23 lbs (10.4kg) 220-240V 50/60Hz, Single Phase, 6.8A

test apparatus for steam jet method ordering information

catalog no.	description	qty
K33900	Existent Gum Evaporation Bath	1
K33850	Series Steam Boiler	1
K33710	Sample Beaker (or 332-002-017)	6
250-000-03F	ASTM 3F Thermometer. Range: +20 to +215°F	2
250-000-03C	ASTM 3C Thermometer. Range: -5 to +400°C	



85 Corporate Drive, Holtsville, New York 11742 1-800-878-9070 (in u.s. only) TEL: +1 631 589 3800 FAX: +1 631 589 3815 Email: sales@koehlerinstrument.com www.koehlerinstrument.com